

Fatty acids are major lipid components, and their content ratios exhibit some features of lipids. In this note, utilizing ADAM (9-anthryldiazomethane) as a fluorescent derivatization reagent, fatty acids are determined by HPLC-fluorescence detection system.

In example 1, the derivatives of representative fatty acids, such as EPA (*cis*-5,8,11,14,17-eicosapentaenoic acid) and DHA (*cis*-4,7,10,13,16,19-docosahexaenoic acid), are successfully separated. However, run time

becomes too long if more hydrophobic fatty acids are analyzed under the condition. Therefore, another condition shown in example 2 was also described in this note.

The derivatization reaction was performed by adding 100 μ L of 0.1 % ADAM solution to 500 μ L of standard solution of fatty acids. After waiting for 1 hour, the reaction solution was injected into the HPLC system.

(K.Suzuki)


Example 1

Conditions


Column :Inertsil ODS-P
(5 μ m, 250 x 3.0 mm I.D.)
Cat.No. 5020-04726

Eluent CH₃CN 0.9 mL/min
Col. Temp. :20 °C
Detection :FL Ex. 365 nm, Em, 412 nm
Inj. Vol. :5 μ L

1.	C10:0	Capric acid	(Decanoic acid)	5 mg/L
2.	C22:5	EPA	(<i>cis</i> -5,8,11,14,17-Eicosapentaenoic acid)	20mg/L
3.	C22:6	DHA	(<i>cis</i> -4,7,10,13,16,19-Docosahexaenoic acid)	5 mg/L
4.	C14:1	Myristoleic acid	(<i>cis</i> -9-Tetradecenoic acid)	5 mg/L
5.	C18:3	Linolenic acid	(<i>cis</i> -9,12,15-Octadecatrienoic acid)	5 mg/L
6.	C12:0	Lauric acid	(Dodecanoic acid)	5 mg/L
7.	C20:4	Arachidonic acid	(<i>cis</i> -5,8,11,14-Eicosatetraenoic acid)	5 mg/L
8.	C16:1	Palmitoleic acid	(<i>cis</i> -9-Hexadecenoic acid)	5 mg/L
9.	C18:2	Linoleic acid	(<i>cis</i> -7,12-Octadecadienoic acid)	5 mg/L
10.	C18:2	Linoleic acid	(<i>cis</i> -7,12-Octadecadienoic acid)	5 mg/L
11.	C16:1	Palmitoleic acid	(<i>trans</i> -9-Hexadecenoic acid)	5 mg/L
12.	C14:0	Myristic acid	(Tetradecanoic acid)	5 mg/L
13.	C18:1	Oleic acid	(<i>cis</i> -9-Octadecenoic acid)	5 mg/L
14.	C18:1	Elaidic acid	(<i>trans</i> -9-Octadecenoic acid)	5 mg/L
15.	C20:1	<i>cis</i> -11-Eicosenoic acid		5 mg/L
16.	C16:0	Palmitic acid	(Hexadecanoic acid)	5 mg/L
		(Deriv.)		

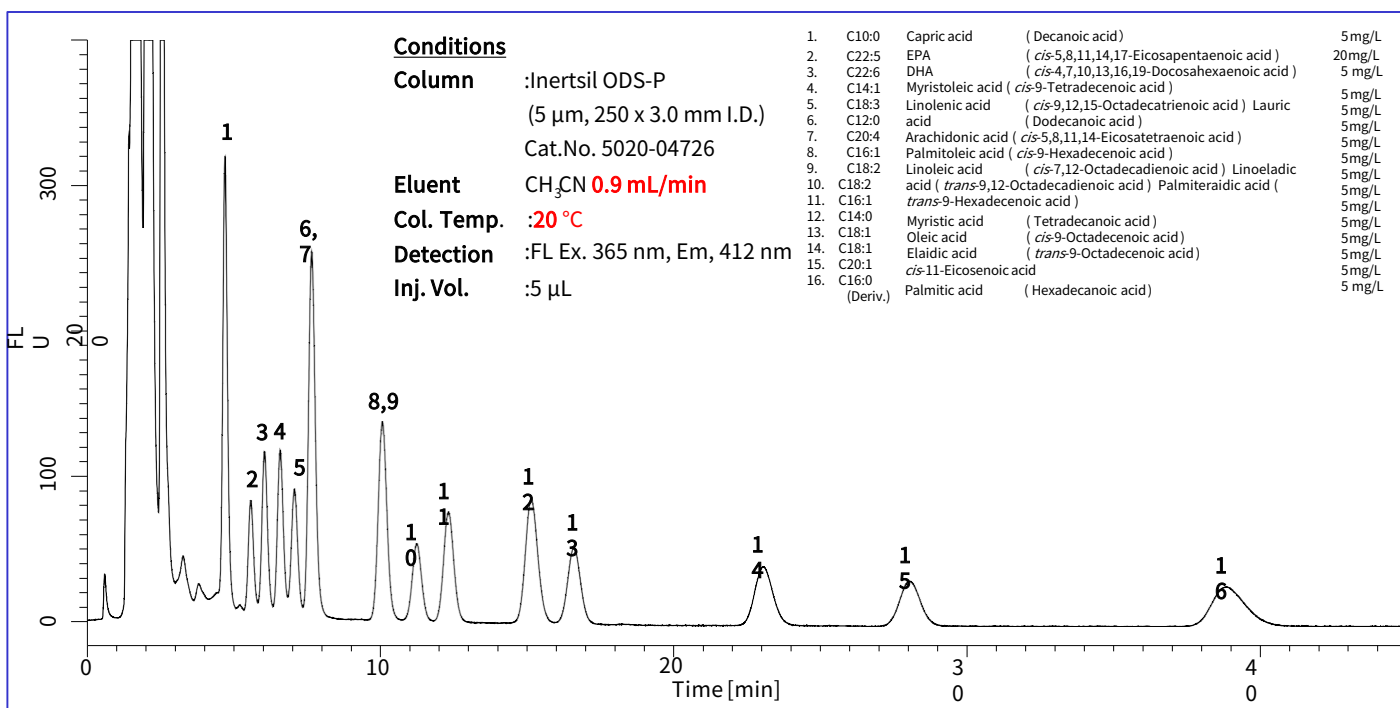

Example 2

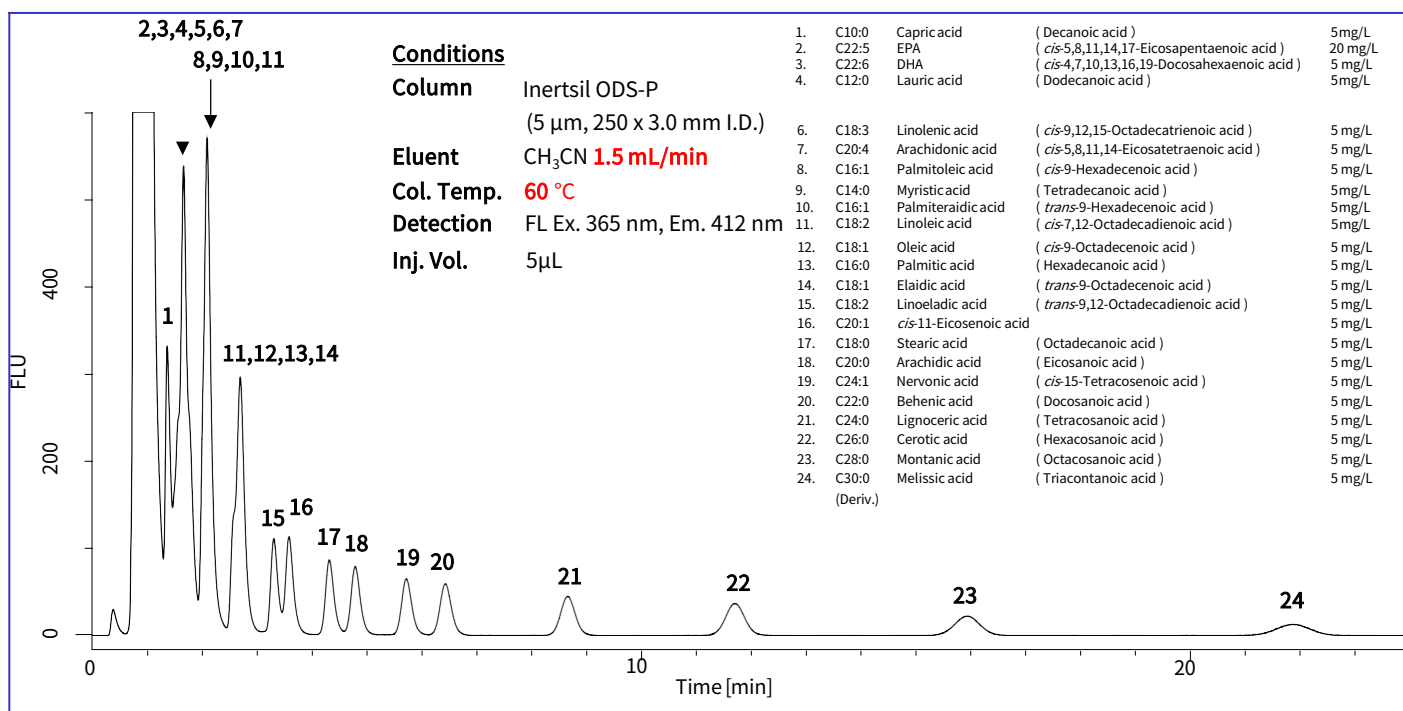

GL Sciences disclaims any and all responsibility for any injury or damage which may be caused by this data directly or indirectly. We reserve the right to amend this information or data at any time and without any prior announcement.

GL Sciences Inc. Japan

22-1 Nishishinjuku 6-chome
Shinjuku-ku, Tokyo
163-1130, Japan

Phone: +81-3-5323-6620
Fax: +81-3-5323-6621
Email: world@glsc.co.jp
Web: www.glsciences.com

GL Sciences Inc. USA

4733 Torrance Blvd. Suite 255
Torrance, CA 90503
USA

Phone: +1-310-265-4424
Fax: +1-310-265-4425
Email: info@glsciencesinc.com
Web: www.glsciencesinc.com

GL Sciences B.V.

Dillenburgstraat 7C
5652AM, Eindhoven
The Netherlands

Phone: +31-40-254-9531
Email: info@glsciences.eu
Web: www.glsciences.eu

GL Sciences (Shanghai) Limited

Tower B, Room 2003
Far East International Plaza
No.317 Xianxia Road, Changning District
Shanghai, China 200051

Phone: +86-21-62782272
Email: contact@glsciences.com.cn
Web: www.glsciences.com.cn



International Distributors

Visit our Website at www.glsciences.com/distributors