


In Japan, a draft of analytical method for linear alkylbenzene sulfonate (LAS) and their salts was released for public comments by Ministry of the Environment. In the draft, enrichment by solid-phase extraction (SPE) and determination using LC/MS/MS are described.

In this note, InertSustain C18 was used as an HPLC column for the analysis. As a result, good reproducibility and linearity of the calibration curve were shown.

(M. Takahashi)

## A Chromatogram Obtained from Standard Solution

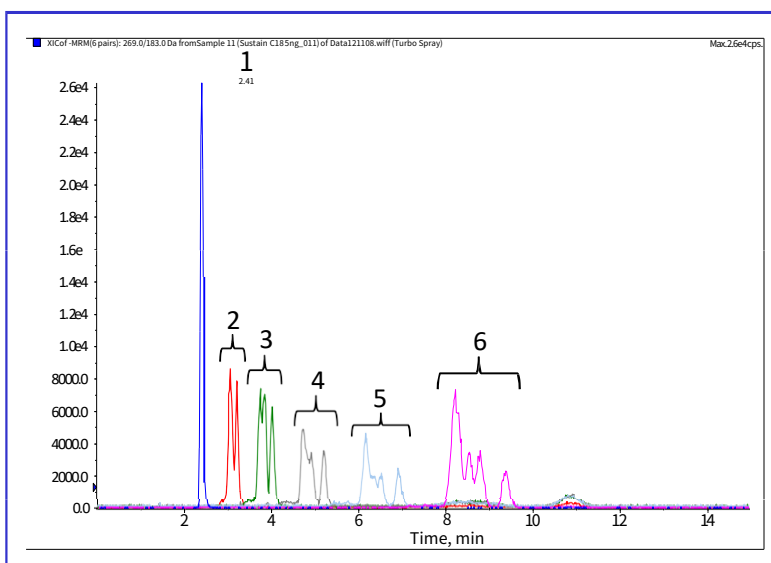

### Conditions

**Column** : InertSustain C18 (3  $\mu\text{m}$ , 150  $\times$  2.1 mm I.D.)

**Eluent** : A)  $\text{CH}_3\text{CN}$

B) 0.1 %  $\text{HCOOH}$ , 50 mM  $\text{HCOONH}_4$  in  $\text{H}_2\text{O}$

A/B = 65/35, v/v

**Flow rate** : 0.2 mL/min

**Col. Temp.** : 40°C

**Detection** : LC/MS/MS

(4000 Q TRAP<sup>®</sup> : ESI, Negative, MRM)

CUR	CAD	IS	TEM	GS1	GS2
10	4	-	600	70	40

**Inj. Vol.** : 5  $\mu\text{L}$

### Analyte:

	Q1/Q3
1. Sodium Octylbenzenesulfonate(C8) (IS)	269/183
2. Sodium Decylbenzenesulfonate(C10)	297/183
3. Sodium Undecylbenzenesulfonate(C11)	311/183
4. Sodium Dodecylbenzenesulfonate(C12)	325/183
5. Sodium Tridecylbenzenesulfonate(C13)	339/183
6. Sodium Tetradecylbenzenesulfonate(C14)	353/183






(in  $\text{CH}_3\text{CN}/\text{H}_2\text{O}$ =65/35 each 5  $\mu\text{g}/\text{L}$ )

### HPLC column


InertSustain C18  
(3  $\mu\text{m}$ , 150  $\times$  2.1 mm I.D.)  
Cat.No. 5020-07415

### Calibration Curves

	Regression equation	Correlation coefficient	RSD, % (50 ng/mL, n=5, Area)
C10-LAS	$y=0.15x-0.0284$	0.9994	1.2
C11-LAS	$y=0.13x+2.31$	0.9988	1.4
C12-LAS	$y=0.117x+1.07$	0.9993	2.7
C13-LAS	$y=0.125x+0.672$	0.9995	2.7
C14-LAS	$y=0.218x+2.56$	0.9975	2.0

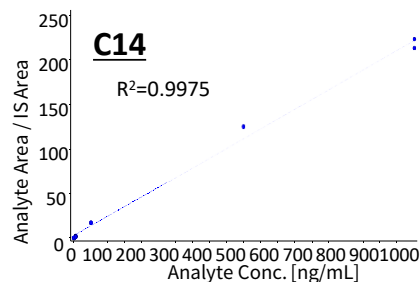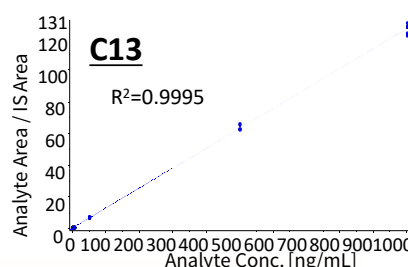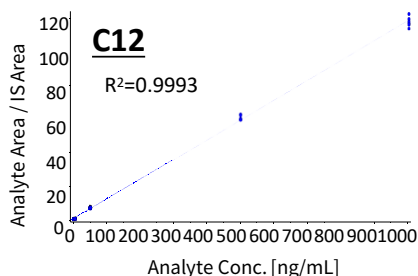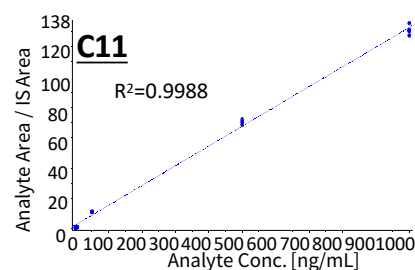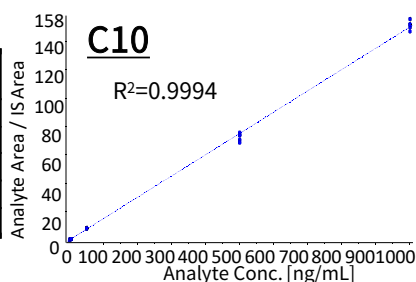

## Another Choice: C8 Column

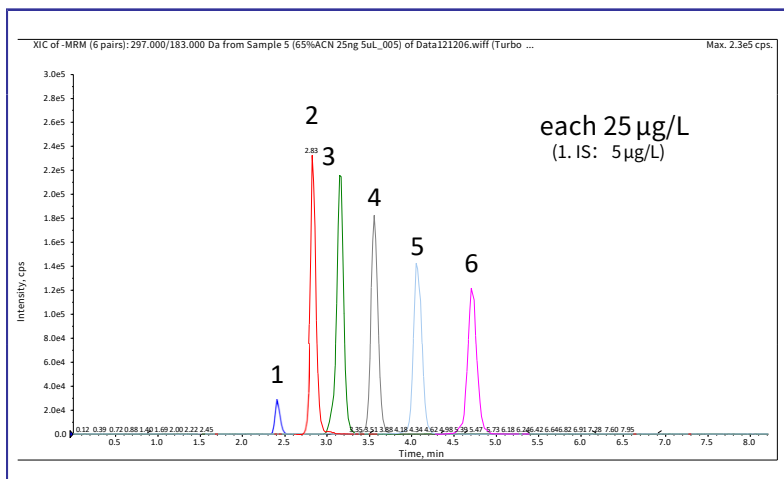


Isomers of each linear alkylbenzene sulfonate are separated to some extent when standard C18 column is used. A chromatogram shown left was obtained by using less retentive C8 column. Each compound was eluted as a single peak because of relatively weak hydrophobic interaction, and peak area can be calculated more easily.

### Conditions

**Column** : Inertsil C8-4 (3 µm, 150×2.1 mm I.D.)  
**Cat.No.** : 5020- 03975  
 Others are the same as described in the previous page.

## Example of Sample Pretreatment using SPE

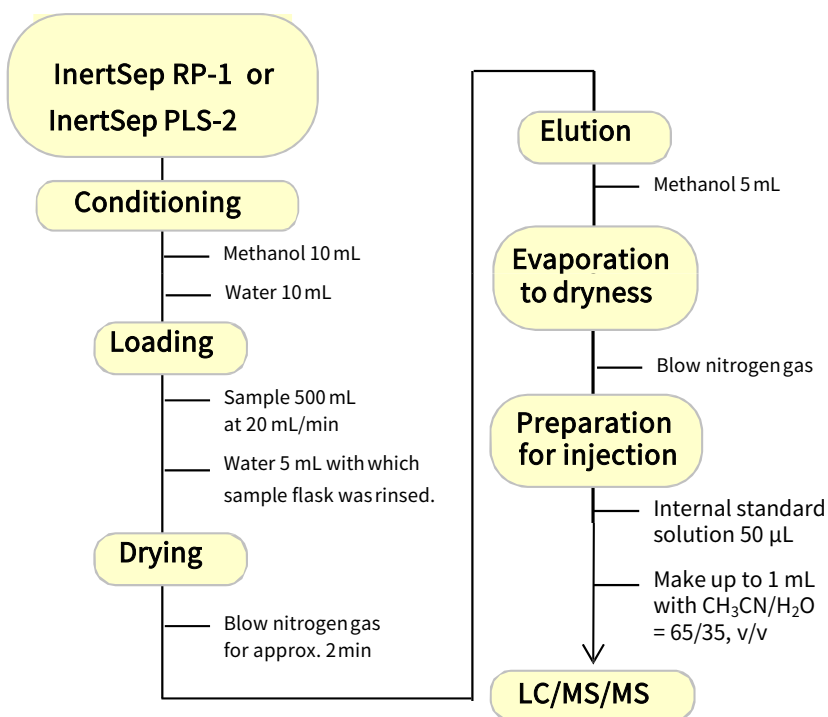

SPE cartridge:  
 InertSep RP-1, InertSep PLS-2



"SlimJ" has top and bottom luer fittings.

- InertSep SlimJ RP-1 230mg 50 pk Cat. No. 5010-65730
- InertSep RP-1 250mg/6mL 30 pk Cat. No. 5010-27000
- InertSep SlimJ PLS-2 265mg 50 pk Cat. No. 5010-65721
- InertSep PLS-2 265mg/6mL 50 pk Cat. No. 5010-27430

GL Sciences disclaims any and all responsibility for any injury or damage which may be caused by this data directly or indirectly. We reserve the right to amend this information or data at any time and without any prior announcement.

**GL Sciences Inc. Japan**  
 22-1 Nishishinjuku 6-chome  
 Shinjuku-ku, Tokyo  
 163-1130, Japan

Phone: +81-3-5323-6620  
 Fax: +81-3-5323-6621  
 Email: [world@glsc.co.jp](mailto:world@glsc.co.jp)  
 Web: [www.glsciences.com](http://www.glsciences.com)

**GL Sciences Inc. USA**  
 4733 Torrance Blvd. Suite 255  
 Torrance, CA 90503  
 USA

Phone: +1-310-265-4424  
 Fax: +1-310-265-4425  
 Email: [info@glsciencesinc.com](mailto:info@glsciencesinc.com)  
 Web: [www.glsciencesinc.com](http://www.glsciencesinc.com)

**GL Sciences B.V.**  
 Dillenburgstraat 7C  
 5652AM, Eindhoven  
 The Netherlands

Phone: +31-40-254-9531  
 Email: [info@glsciences.eu](mailto:info@glsciences.eu)  
 Web: [www.glsciences.eu](http://www.glsciences.eu)

**GL Sciences (Shanghai) Limited**  
 Tower B, Room 2003  
 Far East International Plaza  
 No.317 Xianxia Road, Changning District  
 Shanghai, China 200051

Phone: +86-21-62782272  
 Email: [contact@glsciences.com.cn](mailto:contact@glsciences.com.cn)  
 Web: [www.glsciences.com.cn](http://www.glsciences.com.cn)



**International Distributors**

Visit our Website at [www.glsciences.com/distributors](http://www.glsciences.com/distributors)